

HIV-1 p24 Antigen ELISA

INTENDED USE

The RETRO-TEK HIV-1 p24 Antigen ELISA is supplied for research purposes only. It is not intended for use in the diagnosis or prognosis of disease, or for screening and may not be used as a confirmatory test in diagnostic situations.

The RETRO-TEK HIV-1 p24 Antigen ELISA is an enzyme linked immunoassay used to detect Human Immunodeficiency Virus Type 1 (HIV-1) p24 antigen in research specimens, including cell culture media, as well as human sera and plasma. It can also be used to monitor the purification and biochemical behavior of HIV-1.

Furthermore, the assay may augment or supplant reverse transcriptase measurements traditionally employed to detect the presence of HIV-1. Such enzymatic measurements are not HIV-1 specific.

The RETRO-TEK HIV-1 IC_x/CR_x Kit (ZEPTOMETRIX Catalog #: 0801096) for immune complex dissociation and confirmation of p24 reactivity enhances the detection of p24 antigen in serum or plasma.

Many specimens contain HIV-1 viral loads greater than the range of the standard curve of this assay. The RETRO-TEK HIV-1 p24 Extended Range ELISA (ZEPTOMETRIX catalog # 0801137) will extend the standard curve of this assay to 4 ng/ml.

PRINCIPLE OF THE TEST




Microwells are coated with a monoclonal antibody specific for the p24 *gag* gene product of HIV-1. Viral antigen in the specimen is specifically captured onto the immobilized antibody during specimen incubation. The captured antigen is then reacted with a high-titered human anti-HIV-1 antibody conjugated with biotin. Following a subsequent incubation with Streptavidin-Peroxidase, color develops as the bound enzyme reacts with the substrate. Resultant optical density is proportional to the amount of HIV-1 p24 antigen present in the specimen.

REAGENTS

Materials Supplied:

- **HIV-1 p24 Antibody Coated Microplate for 96 determinations, 1 plate:** 12x8 well strips.
- **HIV-1 p24 Detector Antibody, 1 vial, lyophilized:** Contains biotin-labeled human antibody to HIV-1, human source material non-reactive for antibodies to HIV-1 and non-reactive for HIV-1 p24 antigen, milk, Triton X-100®, Tween 20®, and PBS.
- **Detector Antibody Diluent: 25 ml**
- **HIV-1 p24 Antigen Standard, 0.5 ml:** Contains detergent-disrupted, heat-inactivated viral antigen, goat serum, Triton X-100®, and sodium azide.
- **Lysing Buffer, 5 ml:** Triton X-100® in PBS and 2-chloroacetamide.
- **Streptavidin-Peroxidase, 0.3 ml:** Contains streptavidin conjugated to horseradish peroxidase, PBS, goat serum, and 2-chloroacetamide.
- **Assay Diluent, 100 ml:** Contains goat serum, PBS, Triton X-100®, and 2-chloroacetamide.
- **10X Plate Wash Buffer, 125 ml:** Contains PBS, Tween 20®, and 2-chloroacetamide.
- **Substrate, 0.3 ml:** Contains tetramethylbenzidine (TMB), and dimethyl sulfoxide (DMSO).
- **Substrate Buffer, 50 ml:** Contains citrate/acetate buffer, hydrogen peroxide, and 2-chloroacetamide.
- **Stop Solution, 12 ml:** Proprietary formulation.

PI0801111
Revision: 06
Effective Date: 06/14/2021

REF	Catalog Number		Temperature Limitation
LOT	Batch Code		Expiration Date
RUO	For Research Use Only		Biological Risk

www.ZeptoMetrix.com

ZeptoMetrix LLC • 878 Main Street, Buffalo, NY 14202 USA • Tel (800) 274-5487

This product was manufactured in a facility which has a Quality Management System that is ISO 13485 certified.

PCA# 21-112

ZeptoMetrix®

HIV-1 p24 Antigen ELISA

Catalog Number:
0801111

- **Plate Sealers:** 10
- **Resealable Plastic Bag:** 1

® Triton X-100 is a registered trademark of Union Carbide Chemicals and Plastics Co., Inc. Tween 20 is a registered trademark of Imperial Chemical Industries.

Materials required but not supplied:

- Disposable gloves
- Validated adjustable micropipettes, single and multichannel
- Test tubes and racks for preparing specimen and control dilutions
- Graduated cylinders and assorted beakers
- Validated automatic microplate washer or manual vacuum aspiration equipment
- Validated incubator for 37°C ±1°C
- Validated microplate reader
- Timer
- 1% sodium hypochlorite as disinfectant. May be prepared from household bleach
- Distilled or deionized water

STORAGE

Store all kit reagents at 2° - 8°C. DO NOT FREEZE. When stored properly the kit is stable until the date indicated on the box label.

PRECAUTIONS

- Prior to performing the assay, carefully read all instructions.
- Use universal precautions when handling kit components and test specimens.*
- To avoid cross-contamination, use separate pipette tips for each specimen.
- When testing potentially infectious specimens, adhere to all applicable local, state and federal regulations regarding the disposal of biohazardous materials.
- HIV-1 p24 Antigen Standard contains sodium azide as a preservative. Sodium azide may react with lead or copper pipes to form explosive metal azides. Flush pipes with large quantities of water upon disposal to prevent azide buildup in drains.
- Stop Solution contains hydrochloric acid which may cause severe burns. In case of contact with eyes or skin, rinse immediately with water and seek medical assistance. Wear protective clothing and eye wear.
- Human source material used in the manufacture of the HIV-1 Detector Antibody has been tested and found negative for Hepatitis B surface antigen. The viral lysate used to prepare the HIV-1 p24 Antigen Standard has been inactivated by chemical disruption and heating. Handle these reagents as if capable of transmitting infectious agents.

**MMWR, June 24, 1988, Vol. 37, No. 24, pp. 377-382, 387-38

PREPARATION OF REAGENTS

HIV-1 p24 Detector Antibody: Add 11.0 ml of **Detector Antibody Diluent** to the bottle of **HIV-1 p24 Detector Antibody**. Mix or rock gently for at least 1 hour. Reconstituted reagent can be stored for one week at 2 - 8°C. Alternatively, the reconstituted HIV-1 p24 Detector Antibody can be divided into aliquots and frozen at -20°C until the expiration date of the kit.

HIV-1 p24 Antigen Standard: Prepare a series of **six standards** from the HIV-1 p24 Antigen Standard. Use the dilution scheme in **Table 1**. Any diluted HIV-1 p24 Antigen Standard remaining after the completion of the assay should be disinfected and discarded. Do not save diluted reagent.

PI0801111
Revision: 06
Effective Date: 06/14/2021

REF	Catalog Number	🔑	Temperature Limitation
LOT	Batch Code	🕒	Expiration Date
RUO	For Research Use Only	🚫	Biological Risk

www.ZeptoMetrix.com

ZeptoMetrix LLC • 878 Main Street, Buffalo, NY 14202 USA • Tel (800) 274-5487

This product was manufactured in a facility which has a Quality Management System that is ISO 13485 certified.

PCA# 21-112

HIV-1 p24 Antigen ELISA

Table 1

Preparation of HIV-1 p24 Antigen Standard

Standard Number	Concentration of HIV-1 p24 (pg/ml)	HIV-1 p24 Antigen Standard (µl)	Assay Diluent (µl)
1	125.0	50	950
2	62.5	500 of #1	500
3	31.3	500 of #2	500
4	15.6	500 of #3	500
5	7.8	500 of #4	500
6	0	0	500

Streptavidin-Peroxidase and Substrate: To prepare the Streptavidin-Peroxidase Working Solution and the Substrate Working Solution, use the dilution schemes in **Table 2**. Use Substrate Working Solution within 15 minutes of its preparation. Any diluted Substrate Working Solution and Streptavidin-Peroxidase Working Solution remaining after the completion of the assay must be discarded.

Table 2

Preparation of Streptavidin-Peroxidase Working Solution and Substrate Working Solution

Number of Strips Used	Streptavidin-Peroxidase Working Solution		Substrate Working Solution	
	Streptavidin peroxidase (µl)	Assay Diluent (ml)	Substrate (µl)	Substrate Buffer (ml)
3	40	4.0	40	4.0
6	70	7.0	70	7.0
9	100	10.0	100	10.0
12	120	12.0	120	12.0

Plate Wash Buffer: Dilute **10X Plate Wash Buffer 1:10** in distilled or deionized water prior to use. 1X Plate Wash Buffer may be stored at 2° - 8°C for up to 1 week. Additional bottles of 10X Plate Wash Buffer (ZEPTOMETRIX Catalog #: 0801060) may be ordered.

REF	Catalog Number	🔑	Temperature Limitation
LOT	Batch Code	🕒	Expiration Date
RUO	For Research Use Only	🚫	Biological Risk

HIV-1 p24 Antigen ELISA

TEST PROCEDURE

When assaying serum or plasma samples, sensitivity may be increased by using the RETRO-TEK HIV-1 p24 IC_x/CR_x Kit, (ZEPTOMETRIX Catalog #: 0801096) to dissociate possible immune complexes. Reactive specimens should be confirmed using the IC_x/CR_x Kit confirmatory reagents.

Allow all reagents to reach room temperature before use. Label test tubes to be used for the preparation of standards and specimens. Label each strip on its end tab to identify the strips should they become detached from the plate frame during the assay. If the entire 96 well plate will not be used, remove surplus strips from the plate frame. Place surplus strips and desiccant into the Resealable Plastic Bag, seal and store at 2° - 8°C.

Step 1: Treat specimens in a test tube by pipetting **50 µl** of **Lysing Buffer** into **450 µl specimen** and mix well.

Step 2: Wash each well of the microplate or the selected number of strips 6 times with **300 µl of 1X Plate Wash Buffer** and aspirate. Thoroughly blot by striking inverted microplate or strips on a pad of absorbent towels. Continue striking until no droplets remain in the wells. **Do not allow washed plates to dry completely prior to sample addition. Drying will adversely affect test results.**

Step 3: Leave one well of the microtiter plate empty during the assay. This well is used for a substrate blank. Pipette **200 µl** of **standards #1-6** into duplicate wells.

Step 4: Pipette **200 µl** of each **specimen**, as prepared in **Step 1**, into duplicate wells.

Step 5: Cover microplate with a plate sealer and **incubate** as follows:

- a. Overnight at 37°C ± 1°C, for serum or plasma specimens treated with RETRO-TEK HIV-1 IC_x/CR_x Kit.
- b. Two hours or overnight at 37°C ± 1°C, for untreated serum or plasma, culture fluids or other biological specimens.

Step 6: Aspirate and wash plate as described in **Step 2**.

Step 7: Pipette **100 µl** of **reconstituted HIV-1 p24 Detector Antibody** into each well, except the substrate blank. Cover the microplate with a sealer and **incubate** for 1 hour at 37°C ± 1°C.

Step 8: Aspirate and wash plate as described in **Step 2**.

Step 9: Pipette **100 µl** of the **Streptavidin-Peroxidase Working Solution** into each well except the substrate blank. Cover the microplate with a sealer and **incubate** for 30 minutes at 37°C ± 1°C.

Step 10: Aspirate and wash plate as described in **Step 2**.

Step 11: Pipette **100 µl** of **Substrate Working Solution** into all wells and **incubate** uncovered for **30 minutes at room temperature (18°- 25°C)**. A blue color will develop in wells containing viral antigen.

Step 12: Stop the reaction by pipetting **100 µl of Stop Solution** into each well. A color change from blue to yellow will result.

Step 13: Within **15 minutes**, read the optical density of each well at **450 nm** using a microplate reader.

PI0801111
Revision: 06
Effective Date: 06/14/2021

REF	Catalog Number	⚠	Temperature Limitation
LOT	Batch Code	🕒	Expiration Date
RUO	For Research Use Only	🚫	Biological Risk

www.ZeptoMetrix.com

ZeptoMetrix LLC • 878 Main Street, Buffalo, NY 14202 USA • Tel (800) 274-5487

This product was manufactured in a facility which has a Quality Management System that is ISO 13485 certified.

PCA# 21-112

TEST VALIDITY

Determine the mean optical density values for each standard and specimen. For the test to be valid, it must meet the following criteria:

- The mean optical density of the 0 pg/ml standard and the substrate blank must be less than 0.200.
- The mean optical density of the 62.5 pg/ml standard must be greater than or equal to 0.500

Additional Validity Requirements for the RETRO-TEK HIV-1 IC_x/CR_x Kit

- For immune complex dissociation, the positive control must be above the cut-off for immune complex dissociated specimens and the negative control must be below the cut-off for immune complex dissociated specimens.
- After treatment with the CR_x Neutralization Reagent, the level of the Positive Control must be reduced by at least 50% and the level of the Negative Control must be below the cut-off.

CALCULATION OF RESULTS

Determination of the Cutoff Value:

A predetermined factor is added to the mean of the negative control or negative specimen sample optical value.

- For non-immune complex dissociated serum or plasma samples, and tissue culture samples add 0.030.
- For immune complex dissociated serum or plasma samples add 0.050.

The values above are general guidelines to determine the cutoff value. Alternatively, more accurate results may be obtained by adding at least two standard deviations to the mean optical density from a set of known negative specimen samples to establish the statistical cutoff value.

Samples with absorbance values greater than or equal to the cutoff value are considered positive. Positive results must be confirmed by retesting in duplicate from the same sample source.

Specimens found to be repeatedly reactive for p24 should be confirmed using the CR_x Neutralization Procedure. Refer to Catalog #: 0801096 (RETRO-TEK IC_x/CR_x Kit). For confirmation of reactivity by the RETRO-TEK IC_x/CR_x Kit, calculate the percent reduction of p24 reactivity in controls and specimens using the formula in Figure 1:

Figure 1:

$$\% \text{ Reduction} = 1 - \left[\frac{S_N \text{ (OD)}}{S_C \text{ (OD)}} \right] \times 100$$

Reactivity is confirmed by a reduction of p24 antigen in the specimen by 50% or greater. If reactivity is not confirmed, the specimen is considered indeterminate.

To Quantitate Levels of HIV-1 p24:

Using linear graph paper, plot the concentration of HIV-1 p24 Antigen Standard (pg/ml) on the X-axis versus the mean optical densities for each standard on the Y-axis. Then determine the concentration of HIV-1 p24 antigen in specimens by interpolation from the standard curve. Alternatively, the level of HIV-1 p24 may be calculated by computer using a point-to-point algorithm. Be sure to correct for all dilutions, including the 1 to 1.1 dilution made during the addition of Lysing Buffer.

PI0801111
Revision: 06
Effective Date: 06/14/2021

REF	Catalog Number	🔑	Temperature Limitation
LOT	Batch Code	🕒	Expiration Date
RUO	For Research Use Only	🚫	Biological Risk

www.ZeptoMetrix.com

ZeptoMetrix LLC • 878 Main Street, Buffalo, NY 14202 USA • Tel (800) 274-5487

This product was manufactured in a facility which has a Quality Management System that is ISO 13485 certified.

PCA# 21-112

TYPICAL STANDARD CURVE

HIV-1 Antigen Concentration	Avg. OD at 450nm in overnight assay*
125.00	1.315
62.50	0.840
31.25	0.508
15.60	0.304
7.80	0.184
0.00	0.055

*2 hour assay values may be less than the overnight results. Standard curves may vary as a result of incubation time and temperature, laboratory temperature, etc.




LIMITATIONS OF THE PROCEDURE

Most HIV-1 infected individuals produce antibodies to p24 antigen. The concentration and microspecificity of these antibodies will vary from individual to individual and from bleed to bleed. The observed level of p24 antigen in any specimen containing p24 antibodies may be affected by the following:

- Host antibodies may mask the epitope reactive with capture MAb. Consequently, the optical density may be reduced.
- Host antibodies may mask epitopes that bind to the detector antibody, again reducing the optical density.

After immune complex dissociation and neutralization, host antibodies may recombine with p24 antigen. The recombination of host antibody with antigen is competitive with the capture of p24 by the MAb on the solid phase and thus may mask some of the p24. The rate of recombination is determined by the concentrations of p24 antigen and anti-p24 antibodies, the affinity of the antibodies, as well as temperature and time and is therefore difficult to control. Correlation of results for specimens between repeat test runs and among kits from different manufacturers may be adversely affected by such competitive recombination.

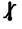PI0801111
Revision: 06
Effective Date: 06/14/2021

REF	Catalog Number		Temperature Limitation
LOT	Batch Code		Expiration Date
RUO	For Research Use Only		Biological Risk


www.ZeptoMetrix.com

ZeptoMetrix LLC • 878 Main Street, Buffalo, NY 14202 USA • Tel (800) 274-5487

This product was manufactured in a facility which has a Quality Management System that is ISO 13485 certified.

PCA# 21-112

PROCEDURAL FLOW CHART

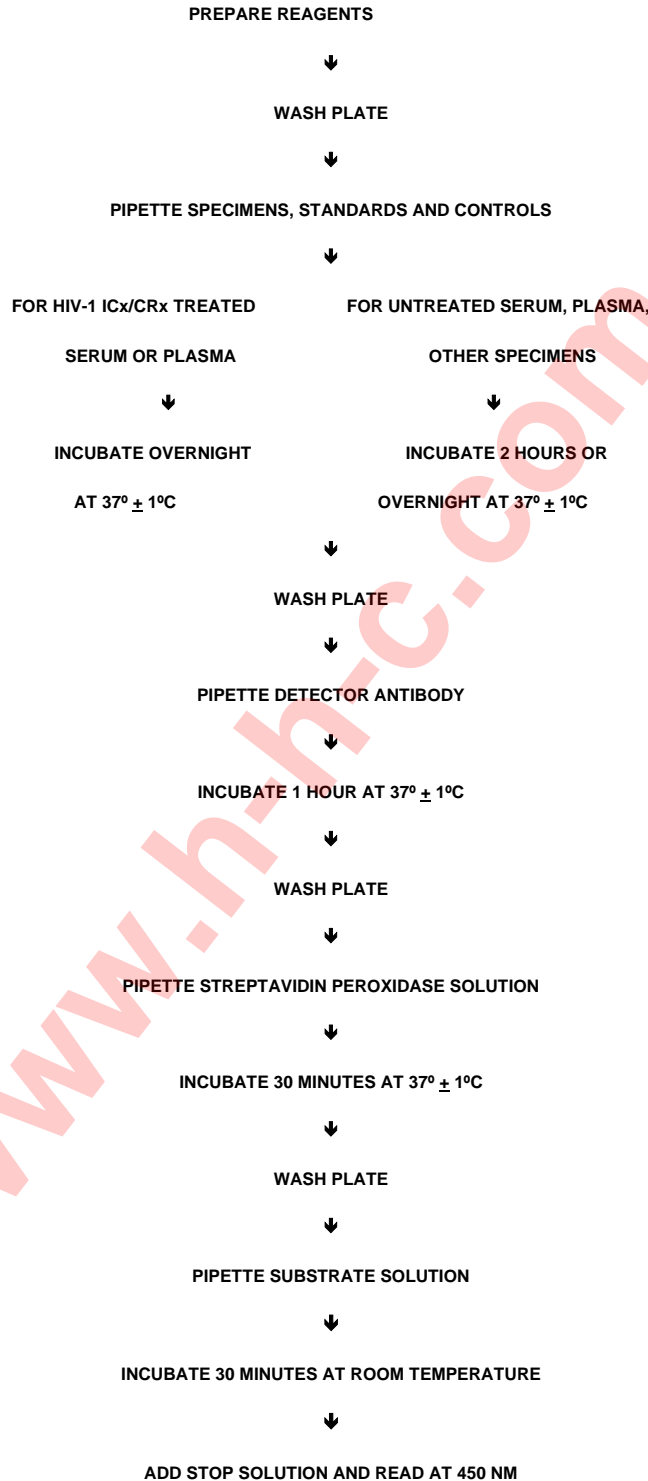

PI0801111
Revision: 06
Effective Date: 06/14/2021

REF	Catalog Number		Temperature Limitation
LOT	Batch Code		Expiration Date
RUO	For Research Use Only		Biological Risk

www.ZeptoMetrix.com

ZeptoMetrix LLC • 878 Main Street, Buffalo, NY 14202 USA • Tel (800) 274-5487

This product was manufactured in a facility which has a Quality Management System that is ISO 13485 certified.

PCA# 21-112